

# Fixation with paraformaldehyde

## Fixation with paraformaldehyde

- Add 41.5ml 2.26% sodium dihydrogen orthophosphate to a suitable heat-proof container
- Add 8.5ml sodium hydroxide 2.52%
- Heat to 60-80°C in a covered container
- Add 2g 'Analar' paraformaldehyde, stirring thoroughly until dissolved
  
- Adding a couple of drops of 1M NaOH can help the PFA dissolve
- Allow to cool and filter to remove particulates. Recheck pH.
- Store at RT and ideally use within 48h.